Pocket Microscope

USER MANUAL







About the Pocket Microscope

The Dartwood Pocket Microscope is compact, portable, and perfect for magnifying tiny objects. It features a powerful magnification range of 60x to 120x, with easy-to-use zoom and focus rings, and a built-in LED light for bright and clear viewing.

In the Box

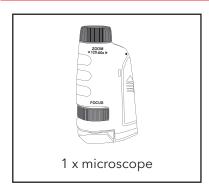

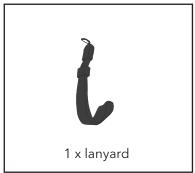


Safety Information

- 1. To avoid permanent eye damage, do not look at the sun or any bright objects with the microscope.
- 2. Remove the battery before storing the microscope for an extended period.
- 3. Inspect all batteries, rechargeable batteries are not recommended. Insert battery in the direction indicated.
- 4. Remove exhausted batteries from the device.
- 5. Small parts on the microscope could be a choking hazard.
- 6. Not suitable for children under 3 years old.
- 7. Children should only use under adult supervision.



Introduction to Parts

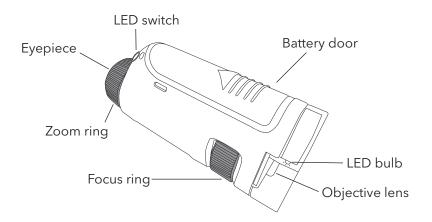


Installing the Battery

1. Slide open the battery door.



2. Install 1 AAA battery (not included) with the polarity shown below.

3. Reattach battery door as shown below.

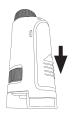


How to Use


1. Place the viewing subject directly below the objective lens.




2. Press the LED button to turn on the LED.

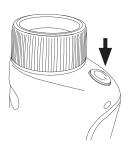


3. Adjust the zoom ring to your preferred magnification.

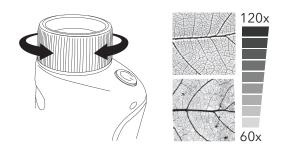

4. Rotate the focus ring until the image is sharp and clear.

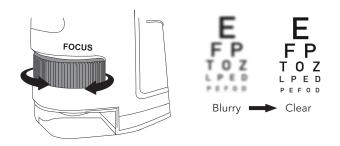

5. Turn off the LED when finished.

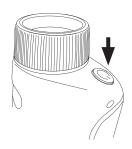

Specifications

Model: DW-KY-A099

Microscope type: 60 - 120x

Focuser: manual

Light source: LED

Materials: plastic Color: white

Battery: 1x AAA (not included) Overall dimension: 3.8 H x 1.8 L x 1.2 W in

Weight: 2.4 oz

Get More Support

Scan the QR code below or email us at contact@wasserstein-home.com







a WASSERSTEIN brance

www.wasserstein-home.com