# Sponge-Stick

## 3M™ Sponge-Stick

English

**REF** see chart.

**Description:** Pre-hydrated and biocide-free 3.8 x 7.6 cm (1.5 x 3 in) sponge in sample bag. Pre-hydrated with specified agent. Various bag sizes. With or without gloves.

**Applications:** For Laboratory Use Only. Material Safety Data Sheet available.

### Product identity:

| Product identity:                                     | Volume | REF        | Volume of Bag  | Gloves |  |
|---|--------|------------|----------------|--------|---|
| 3M™ Sponge-Stick                                      | 0 mL   | SSL100     | 0.89 L (30 oz) | 0      | 2-30°C  |
| 3M™ Sponge-Stick with<br>Buffered Peptone Water Broth | 10 mL  | SSL10BPW   | 0.89 L (30 oz) | 0      | 2-30°C  |
| 3M™ Sponge-Stick -<br>Lethen Broth - 2 Gloves         | 10 mL  | SSL10LET2G | 0.89 L (30 oz) | 2      | 2-8°C   |
| 3M™ Sponge-Stick -<br>Neutralizing Buffer             | 10 mL  | SSL10NB    | 0.89 L (30 oz) | 0      | 2-30°C  |
| 3M™ Sponge-Stick -<br>Neutralizing Buffer - 2 Gloves  | 10 mL  | SSL10NB2G  | 0.89 L (30 oz) | 2      | 2-30°C  |


### Instructions for Use

- 1) Shake stick to end of bag.
- 2) Tear bag open.
- 3) Squeeze bag to open. Aseptically grasp the stick above the thumb-stop line to remove the sponge.
- 4) Aseptically swab across the entire sampling surface.
- 5) Turn sponge over. Change direction 90°. Aseptically swab the same sampling surface.
- 6) Aseptically place sponge into bag up to the thumb-stop.
- 7) Hold sponge in place inside bag. Bend stick to break. Allow sponge to drop in bag. Discard stick.
- 8) Fold bag to close.
- 9) Fold ends of blue wire inward.

**Disposal:** Follow established laboratory practices.

Storage Symbol: See chart.

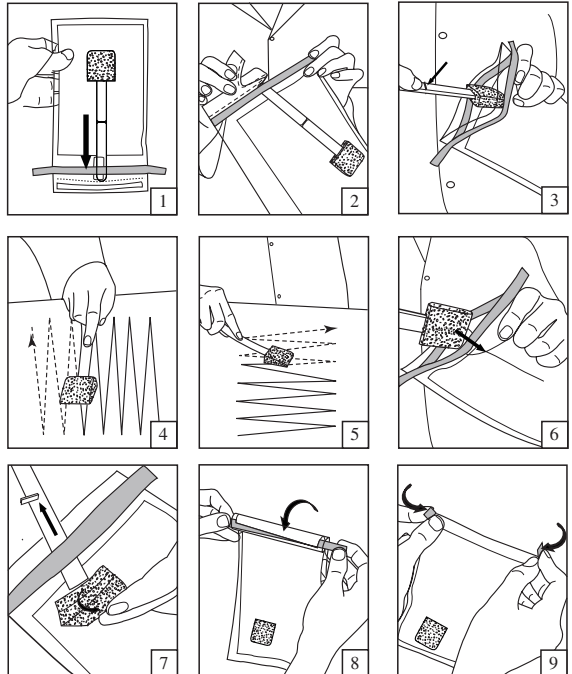## 3M

St. Paul, MN, 55144-1000 USA  
www.3M.com/microbiology

3M is a trademark of 3M and used under license in Canada. ©3M 2008.  
All rights reserved.  
Made in U.S.A.

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©3M 2008. All Rights Reserved.