

Avalon International Corp.

317 Lockwood Avenue,
Phone: 847-446-9986
Web: www.bugcheck.us

Northfield, IL 60093
Fax: 847-446-9946
Email: info@bugcheck.us

BUG CHECK® FC BACTERIA CONVERSION TABLE

Interpretation Chart Value	0	1	2	3	4	5
LIQUID SAMPLE						
Bacterial Count/ml-Undiluted	Less than 100	100	1,000	10,000	100,000	1 million
Bacterial Count/ml-1:10 Dilution	Less than 1,000	1,000	10,000	100,000	1 million	10 million
Bacterial Count/ml-1:100 Dilution	Less than 10,000	10,000	100,000	1 million	10 million	100 million
Bacterial Count/ml-1:1000 Dilution	Less than 100,000	100,000	1 million	10 million	100 million	1 billion or more
SURFACE SAMPLE						
Bacterial Count/in2	Less than 10	10	100	1,000	10,000	100,000

BUG CHECK® FC MOLD OR YEAST CONVERSION TABLE

Interpretation Chart Value	1	2	3	4	5
LIQUID SAMPLE					
Mold or Yeast Count/ml-Undiluted	Less than 10	10	100	1,000	10,000 or more
Mold or Yeast Count/ml-1:10 Dilution	Less than 100	100	1,000	10,000	100,000 or more
Mold or Yeast Count/ml-1:100 Dilution	Less than 1,000	1,000	10,000	100,000	1 million or more
SURFACE SAMPLE					
Mold or Yeast Count/in2	Less than 1	1-3	10	100	1,000 or more

BUG CHECK® FC

Test System for Counting Aerobic
Bacteria and Fungi in Fuel



Contains all supplies necessary to perform
40 tests: 20 bacteria tests + 20 fungi tests

Read all directions before using this product.
Store in refrigerator for maximum shelf life.