



Biocrede Incorporated

Tel: 1-734-678-6956
Email: info@biocrede.com
Website: www.biocrede.com

TaqExpress DNA Polymerase

Store at -20°C

Cat. No.	Description	Quantity
PCR-0082	TaqExpress DNA Polymerase (1000U)	200 µl
PCR-0081	TaqExpress DNA Polymerase (250U)	50 µl

Product Description

TaqExpress DNA polymerase is an engineered version of Taq DNA polymerase developed to achieve rapid PCR. It catalyzes the 5'-3' synthesis of DNA. This enzyme possesses 5'-3' exonuclease activity and moderate 3'-5' proofreading exonuclease activity. The extension speed is about 6 kb/min, which is 6 times faster than the regular Taq DNA polymerase. Template-independent "A" can be attached at the 3' end of the PCR product which can then be cloned into a TA cloning vector.

Product Components	250U	1000U
TaqExpress DNA Polymerase (5 U/µl)	50 µl	200 µl
5X PCR Buffer, with Mg ²⁺	1 ml	4 ml
25 mM MgSO ₄	1 ml	1 ml

Unit Definition

One unit of the enzyme catalyzes the incorporation of 10 nmol of deoxyribonucleotides into a polynucleotide fraction in 30 mins at 74°C.

Storage Buffer Components

50 mM Tris-HCl (pH 8.0), 100 mM NaCl, 0.1 mM EDTA, 5 mM DTT, 50 % glycerol and 1.0 % Triton[®]X-100.

Shipping and Storage

Upon arrival, TaqExpress DNA Polymerase should be stored at -20°C. Avoid repeated freeze-thaw cycles of all TaqExpress components to retain maximum performance. All TaqExpress components are stable for 1 year from the date of shipping if stored and handled properly.

Protocol

The following basic protocol serves as a general guideline and starting point for any

PCR amplification. Optimal reaction conditions (incubation times, temperatures, concentration of Taq DNA Polymerase, primers, MgSO₄ and template DNA) may vary and need to be optimized for each specific PCR.

All PCR experiments should be assembled in a nuclease-free environment. In

addition,

DNA sample preparation, reaction set-up and subsequent reaction(s) should be performed in separate areas to avoid cross contamination.

A negative control reaction (omitting template DNA) should always be performed in tandem with sample PCR to confirm the absence of DNA contamination.

1. Add the following components to a sterile 0.2 ml PCR tube sitting on ice.

Components	Volume	Final Concentration
Template DNA	~100 ng	~2 ng/µl
Forward primer (10 µM)	1 - 2.5 µl	200 - 500 nM
Reverse primer (10µM)	1 - 2.5 µl	200 - 500 nM
5X PCR Buffer, with Mg ²⁺	10 µl	1X
25 mM MgSO ₄ (optional)*	0 - 3 µl	1.5 - 3 mM
dNTP Mix (10 mM)	1 µl	200 µM
TaqExpress DNA Polymerase (5 U/µl)	0.5 - 1 µl	2.5 - 5 U
Nuclease-free H ₂ O	up to 50 µl	-

Notes: Optimal Mg²⁺ concentration is specific to each DNA template-primer set and can only be determined experimentally. We recommend preparing a master mix for multiple reactions to minimize reagent loss and enable accurate pipetting.

2. Mix contents of tube and centrifuge briefly.
3. Incubate tube in a thermal cycler at 94°C for 3 mins to completely denature the template.
4. Perform 30 - 35 cycles of PCR amplification as follows:
Denature: 94°C for 5 sec
Anneal: 45 - 72°C for 15 sec
Extend: 72°C for 10 sec/1 kb template
5. Incubate for an additional 5 mins at 72°C and maintain the reaction at 4°C. The samples can be stored at -20°C until use.

Analyze the amplification products by agarose gel electrophoresis and visualize by ethidium bromide or Non-Toxic Green (Cat No. PCR-7011-NT) staining. Use appropriate molecular weight standards.

*For laboratory research only. Not for clinical applications.
For technical questions, please email us at technical@biocrede.com or
visit our website at www.biocrede.com*