



Vacc-Safe® RANGE

VS250P

VS400P

VS600P

VS1000P

READ THIS BEFORE PLUGGING IN UNIT

Congratulations on choosing a Vacc-Safe® Premium Medical Refrigerator. This set of instructions will show you how to properly install and set up your Vacc-Safe® unit to ensure optimal performance. Failure to follow these instructions may result in a **Service Charge** if a technician needs to attend.

1. Remove all packaging and wrap from the Unit
2. Check for any damage to the unit. Report any damage immediately
3. **INSIDE THE UNIT:**



REAR SPACER



2 x KEYS



SHELVES



OR




SHELF CLIPS



QR CODE

(The VS250 has a metallic U-shaped bracket)

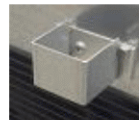


4. Fit the REAR SPACER provided by using the 2 x screws provided (requires Philips screwdriver *not supplied*). This prevents the Unit being pushed hard against a rear wall and allows the Condensing Unit adequate ventilation.

Failure to fit the REAR SPACER will **VOID the WARRANTY.**


5. Keep the unit on a firm, flat and even surface. Do not place the unit in direct sunlight or near heat sources. Ensure the room has adequate ventilation.
6. Once the Unit has been positioned in its location, the front Castors have a BRAKE. Lock the brake on both sides to prevent the Unit from moving when pulling on the handle to open the door.



7. **DOOR(s):** The door(s) on the Vacc-Safe Refrigerator is Lockable. Insert the key serrated side 'down' to Lock / Unlock the Door. Keep one key in a safe place as replacement keys incur a fee.
8. **SHELVES:** Position the shelf clips at even heights and regular intervals suitable for the intended contents of the Unit and place the shelves once finished. The MAX LOADING per shelf is 15kg.
DO NOT STORE PRODUCT DIRECTLY ON THE BASE OF THE UNIT
 - a) Quantity of Shelves per unit
 - i) VS250P x 5
 - ii) VS400P x 7
 - iii) VS600P x 6
 - iv) VS1000P x 12
9. **DO NOT** place items on the shelves in such a way that the entire surface area is covered. This will lead to restricted airflow and reduce the efficiency and effectiveness of the Unit. A recommendation of **¾ maximum usage of the width, height and depth** will allow the unit to function properly.
10. **START-UP:** Once in position and levelled, plug in the unit to a normal GPO (General Power Outlet).



If the Display reads Off, 'press and hold' the power button  on the Control Fascia for 5 seconds; the unit will start within a few seconds and the Cooling Cycle will begin.

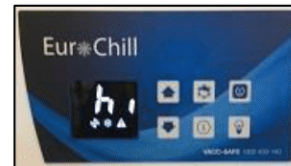11. **On initial start-up**, depending on the ambient temperature, the Unit may temporarily go into 'Hi-Temp' Alarm Mode.
 - To silence the Alarm, press ANY KEY.



12. To turn the internal light on or off use the switch 
13. Once the Unit has reached approximately 5°C you may begin to store items inside. Ideally, they will already be at the desired temperature. Once the unit is in range perform a reset of the high and low temperature readings. The Unit has a QR CODE next to the front control panel which will direct you to a YouTube video instruction guide.
 - a) **NOTE:** Loading the unit with room temperature product will cause the ambient temperature in the fridge to exceed 8°C and cause an alarm. It will also put excessive strain on the compressor and shorten the life of the Unit.
14. Initialise and start your DATA LOGGER as per the instructions included with the unit.
15. The INSTRUCTION MANUAL for this unit can be located by scanning the QR CODE  located on the Model Number label on the inside of the Unit.

If you have any Questions, please call our Help Line – 1300 459 140