
# LIGHTTELLS Corporation Limited Coffee Caffeine Chlorogenic Acid Analyzer

**CA-700** 

**User Manual** 



## Operation Instruction

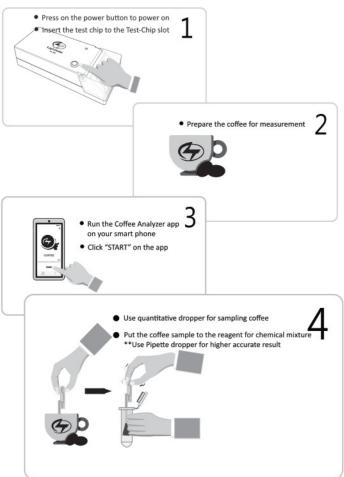

### Product Introduction

The CA-700 portable Coffee Caffeine / Chlorogenic Acid (CGA) Analyzer is an accurate instrument for measuring the caffeine and CGA of coffee with its advanced mobility, accuracy, easy-to-operate feature, you can get the Caffeine and CGA value displayed on your smart phone app within 1 minute.

### Package Contents

- CA-700 x1
- 5V Power Adapter x 1
- USB type C charging cable
- Test Chip package x 10
  - Test Chip x 1
  - Reagent x 1
  - Dropper x 1 (for mixing reagent)
- User Manual with Limited Warranty Page x 1

### Major Features

- One-click quick measurement
- Premix reagent
- Result in 1 minute
- Bluetooth connec@on for iOS/Android app

### ninute

# • Shake the mixture for at least 10 seconds • Use another dropper for the mixture • Drop the mixture on the reacting area of the Test Chip • Make sure the reacting area is fully covered with the mixture

Waiting the result display on the app

Chlorogenic: CGA mg/dl

Caffeine: CAF mg/dl

### Specification

Reading Range	Caffeine 10 – 500 mg/dL
	CGA
Sampling Size	0.1ml
Dropper tolera	ance * Pipette +/ - 6% (25 degree C)
Dropper tolera	ance * Dropper +/- 12% (25 degree C)
Working Hour	2 – 3 conĢnuous working hour per full ch
Weight	370g
Operating temperature	15~35°C
Storage temperature	5~45°C
Dimension	L: 202mm
	W: 78mm
	H: 40mm
Power adapter	110-240V, 50/60Hz
Charging input	5V, 0.5A, USB Type C

### App installation

- iOS: Search "Coffee Analyzer" on App Store
- Android: Search "Coffee Analyzer" on Google Play



### Classification of hazardous info.



**Substance: Coffee detection reaction solution** 

Hazardous ingredients: Sulfuric acid, oxalic acid, citric acid

Hazard warning message: Warning

Health hazard: Corrosive/irritating skin substances

### Hazard prevention:

Please operate the product in a well-ventilated place. It is strictly Forbidden from eating or drinking during the testing process, If the reagent contacts the eyes or skin, promptly wash (irrigate) the eyes with large amounts of water or seek medical attention as soon as possible.

### Manufacturer:


(1) Company: Lighttells Corp.. Ltd.


(2) Address: 5F, No. 27, Xienzheng 3rd St., Zhubei city,

Hsinchu county, Taiwan, R.O.C.

(3) TEL: +886-3-555-1389


### CA-700 Hardware

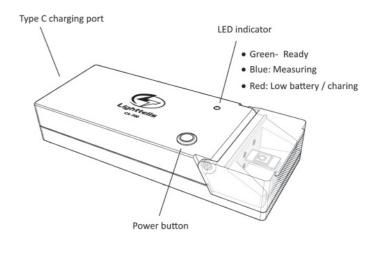

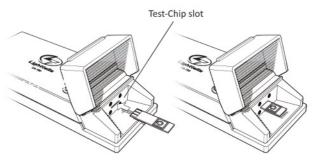

Side up the test-chip, plug into the tes@ng slot

### Package Contents

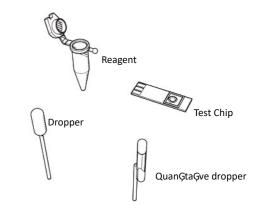

### Notice

- The tesGng pack is for Single-Use only, reuse IS NOT RECOMMENDED.
- For stable accurate result, it is not recommended charging while tesĢng. Please stay the CA-700 on a stable plane.

### **Warranty Information**

This Limited Warranty applies to LIGHTTELLS product model name CA-700 purchased from an authorized LIGHTTELLS dealer by the original purchaser for normal use and not for resale. LIGHTTELLS warrants that a CA-700 product is free from defects in materials and workmanship, with the exceptions stated below.

This limited warranty lasts for 1 year from date of purchase, depending on when and where it was originally purchased. A valid proof of purchase may be required to prove eligibility. If you do not have a valid proof of purchase, the limited warranty period will be measured from the date of sale from LIGHTTELLS to the authorized distributor.

The limited warranty does not cover damage resulting from improper or unreasonable use or maintenance, misuse, accident, modification or alteration to hardware or software, tampering, unsuitable physical or operating environment beyond product specifications, or failure caused by a product for which LIGHTTELLS is not responsible. There is no warranty of uninterrupted or error-free operation. There is no warranty for product with removed or altered identification labels. LIGHTTELLS DOES NOT PROVIDE ANY OTHER WARRANTIES OF ANY KIND. INCLUDING, BUT NOT LIMITED TO, THE IMPLIED WARRANTIES OR CONDITIONS OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. SOME JURISDICTIONS DO NOT ALLOW THE LIMITATION OF IMPLIED WARRANTIES, SO THIS LIMITATION MAY NOT APPLY TO YOU, LIGHTTELLS is not responsible for returning to your product which is not covered by this limited warranty.

### Limited Warranty Card

Model Name: CA-700 Coffee Caffeine Chlorogenic Acid Analyzer

Serial Number: YY

Date of Purchase: MM DD

Customer Name:

Customer Address:

Authorized Dealer Name / Stamp:

