

20X ELECTROPHAST BUFFER + CONCENTRATE

950ml +  = 1L
of Di Water of ElectroPhast 1X Running Buffer


IBI's 20X Super Fast Agarose gel running buffer was developed to run agarose gels 20–30% faster than standard running buffers, while also providing sharper and better separated bands!

Actual Size

Product Number	Size
IB70190	50ml (20X Solution)
IB70191	1L (20X Solution)


Key Features:

- Use IBI ElectroPhast for a running buffer and for making agarose gels
- Run higher voltages and faster gels: 500V–600V for 15min.
- Get clearer bands and better separation
- Run downstream procedures as always. Does not affect gel extraction or subcloning.
- A great value! As little as \$1.55 per/Liter for running buffer!

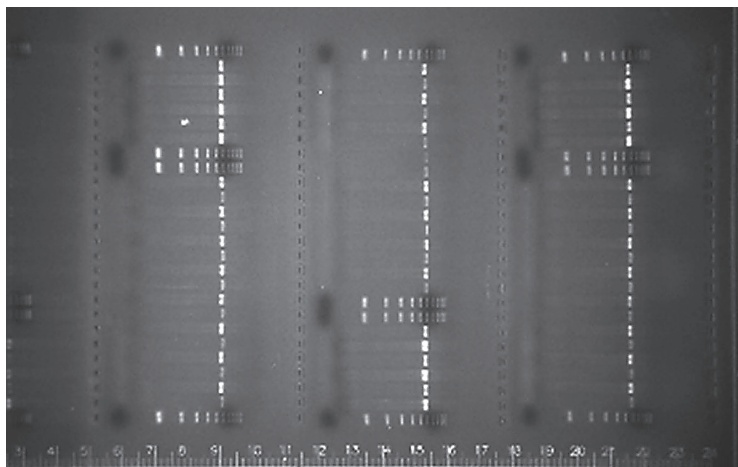

IBI ELECTROPHAST BUFFER AT 600V FOR 15MIN.

- Note** the bands are straight
- Note** the minimal streaking
- Note** the good separation at High MW

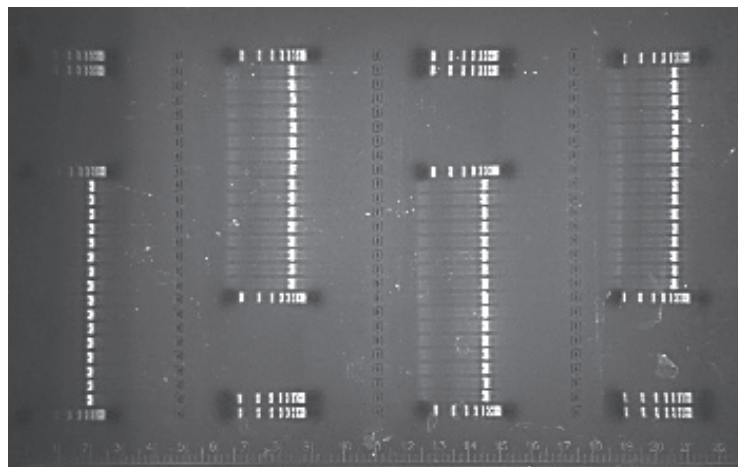

10X TBE GEL RUN AT 600V FOR 15MIN.

- Note** the band cupping & overloading
- Note** the streaking
- Note** the separation at High MW



SCIENTIFIC

800-253-4942 • IBISCI.com ISO 9001 Certified

Building Quality Scientific.