



IBI SCIENTIFIC

9861 Kapp Court Peosta, IA 52068

Tele: (563) 690-0484 Toll Free: 1-800-253-4942

Fax: (563) 690-0490 E-Mail: info@ibisci.com

Web: www.ibisci.com

ISO 9001:2008 Certified

ACCAA0002 – The Belly Dancer Nesting Platform Installation

Important! – Prior to starting project, ensure unit is unplugged from electrical source.

- 1.) To install The Belly Dancer Nesting Platform, you must first remove the existing Belly Dancer Platform. To do this remove the (3) phillips drive screws located at the center of the platform. Next, locate the four flush-mounted allen head screws attaching the platform to the flexible legs. Use an allen wrench to remove these (4) screws. The existing Belly Dancer Platform is now free, and can be removed from the unit.
- 2.) Now, locate the new nesting platform and hardware to prepare for installation. Place the nesting platform onto the top of the Belly Dancer and align the three holes in the center of the platform with the holes located at the top of the tilt assembly below. Insert the (3) phillips head screws and tighten. It works best to start all three screws first, and then tighten them after all screws are mated to the tilt assembly. Next, align the fitting at the top of the flexible leg with the hole at each corner of the nesting platform. Insert the allen head screw and thread into the fitting, as you tighten this screw you will feel the fitting draw up into the acrylic. This is what secures the two parts together, continue to tighten until fitting is fully seated into acrylic and screw is tight.
- 3.) Finally, place new white nesting platform mat onto platform.
- 4.) Reconnect unit's power cord to input entry module and standard power source. Turn control dial on front of unit and ensure unit will now function properly.

