



EJ-07 / EJ-08 Underhook

By attaching the underhook to the bottom of the balance, large objects that are difficult to load on the weighing pan can be weighed, and the density (specific gravity) of objects may be measured. Refer to the "9-6. EJ-13 Density Determination Kit" for the information about the density measurement.

- ❑ ***EJ-07 is for use with the EJ-3000 / EJ-4100 / EJ-6100.***
- ❑ ***EJ-08 is for use with the EJ-1500 / EJ-2000.***
- ❑ ***The calibration with a weight being hung on the hook is required for an accurate weighing.***

EJ-07 / EJ-08 Installation

Open the cap on the bottom of the balance, and screw the underhook into the mounting hole.

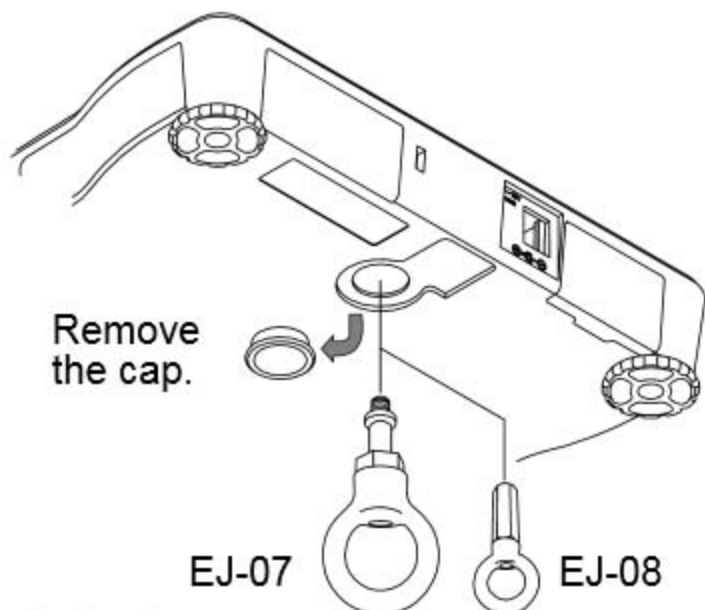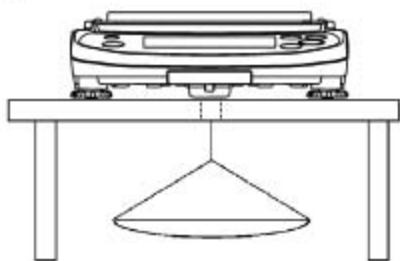

Caution

- ❑ ***Do not apply excessive force to the underhook.***
- ❑ ***When not in use, remove the underhook and attach the cap to prevent dust from getting into the balance.***