

**Serum-Free Recombinant RocketFuel™  
(SF-4ZR-500-R)**

**Product Description:** RocketFuel-R™ Supplement contains recombinant human growth factors and has been designed to provide defined growth supplements in the serum-free environment. The 8mL vial activates 500mL Cell Systems Serum-Free Medium, SF-4Z0-500-R. RocketFuel-R™ is the broad-spectrum mitogen growth supplement designed for use with Serum-Free Medium. RocketFuel-R™ is also available individually as a stand-alone supplement to improve performance of your serum-free medium.

Cell Systems media and reagents are sterile, made with WFI and all components are cGMP and ISO Compliant.

**Components:** This 8mL vial supplements 500mL Serum Free Medium. RocketFuel-R™ is included in Cell Systems Cell Culture Serum Free Medium Kit, SF-4Z0-500-R.

**Storage:** Can be frozen at -20°C. Once opened, shelf life is 30 days at +2-8°C. If entire unit will not be used within 30 days, aliquot and freeze.

**Product Use:** SF-4ZR-500-R is for laboratory research only. It is not approved for human or animal use, or for application in *in vitro* diagnostic procedures.

**Shipping:** All medium and reagents shipped at ambient temperature.

**Instructions for Use:** Make sure all work surfaces are disinfected. Spray the outside of your gloves prior to work under the hood.

- Spray the outside of all tools/instruments (bottles, tubes, racks, etc.) with ethanol solution before bringing them under the hood. Use 8mL RocketFuel-R™ to supplement 500mL Serum Free cell culture medium, or our SF-4Z0-500-R Serum Free Medium.
- Open caps on Serum Free medium and RocketFuel-R™, place caps on gauze pads saturated in alcohol.
- Pour or pipette 8mL RocketFuel-R™ into 500mL Serum Free Medium, SF-4Z0-500-R.