Caframo

Petite Digital Stirrer User Manual

Online Warranty Registration

To register the warranty for your Caframo stirrer Online, please use the simple registration form on our website.

www.CaframoLabSolutions.com

Mail or Fax Warranty Registration

To register the warranty for your Caframo stirrer via Mail or Fax, please fully complete the reverse form. Once completed, detach this page from the manual and return to us using one of the following methods:

Mail

Attn: Caframo Lab Solutions - Warranty Validation Department
501273 Grey Road 1
Georgian Bluffs ON NOH 2T0
Canada

Fax

800-209-6786

Caframo Warranty Registration

Company Name _____ Contact Name _____

Address _____ City _____ State/Prov _____ Country _____ Zip/Postal Code _____

Telephone _____ Fax _____ E-mail _____

Name of Dealer that supplied your Caframo Product _____ Date of Purchase _____

Model # _____ Serial # _____

Please note the correct responses and return to us via the instructions on the reverse.

A. Frequency of use:

___ Continuous Duty (1)

___ Approx 8 hrs/day (2)

___ Approx. 4 hrs/day (3)

___ Approx. 1 hr/day (4)

___ Infrequent Use (5)

C. Volumes mixed:

___ 0-1 litre (1)

___ 1-2 litres (2)

D. Average speed range used:

___ 1-1000 RPM (1)

___ 1000-2000 RPM (2)

F. Do you use other Caframo Products?

___ Yes ___ No

G. Do you use similar products from other manufacturers?

___ Yes ___ No

If yes, who is the manufacturer? _____

B. Viscosity of product:

___ Low Viscosity

(1CPS - ie. Water) (1)

___ Med. Viscosity

(1000CPS - ie. Paint, Oil) (2)

E. Why did you purchase this product?

___ Reputation (1)

___ Pricing (2)

___ Capabilities (3)

___ Other (4)

I. What do you require in your facility that you are having a difficult time obtaining?

Language Index

English.....	1
French.....	9
Spanish.....	17
German	25

Contents

Language Index	1
Packing List	1
Safety Instructions	2
General Instructions	3
Approvals	3
Description	4
General Operating Instructions	4
Technical Data	5
Environmental Operating Conditions	5
Installation in the Laboratory	6
Cleaning and Service	6
Warranty and Product Liability.....	7
Accessories.....	7

Packing List

Compare the contents of the shipment (BDC250 or BDC250U1) to the items on the following list to ensure all parts are received with the stirrer. Do not discard the container and packing materials until all parts are accounted for. It is recommended that the original packing materials be kept for transportation and storage.


- Petite digital stirrer *
- Power supply *
- Instruction manual *
- Allen key, (2) collet set screws *
- Stand support rod and lock nut
- Stand base
- (2) Impeller blades
- (2) Impeller lock screws
- Impeller shaft

* Included with BDC250MU1 (motor only)

Universal Power Supply

Included with BDC250U1, BDC250MU1.

- Interchangeable plugs for:
 - a. North America/Mexico/Japan
 - b. Europe
 - c. Australia/China
 - d. UK/Singapore
- Measures H 74, W 43.5, D 34mm
- RoHS compliant & CE marked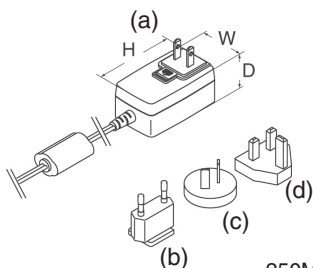

Safety Instructions

- ⚠ 1. Each operator should read all instruction before using this device.
- ⚠ 2. The stirrer must be securely fixed to a stable support such as the provided stand. If other stands are used, the unit must be secured such that it will not fall if tipped.
- ⚠ 3. **CAUTION:** This is not an explosion proof stirrer. Do not use in a highly flammable or explosive environment.
- ⚠ 4. Spinning impellers can cause severe personal injuries. Operators must use extreme care and good judgment when mixing at any speed.
- ⚠ 5. All mixing paddles and impellers must be in good condition with straight shafts. If the stirrer vibrates at high speeds, check the paddle shaft for damage, and repair or replace if needed.
- ⚠ 6. Extreme care must be taken when mixing chemicals to ensure that no chemicals are splashed outside the mixing vessel. Care must be taken when changing to faster mixing speeds.
- ⚠ 7. Ensure that the mixing impeller does not contact the containment vessel.
- ⚠ 8. Do not operate while standing in water. Keep the unit dry and do not immerse any part, except the mixing paddle into any liquids. Protect unit and operator from splashing.
- ⚠ 9. Ensure that no loose clothing, jewellery, or hair can become entangled in any rotating parts. A fast spinning chuck can cause injury to operator.
- ⚠ 10. Power can be interrupted to the stirrer by pressing the power button. In an emergency, disconnect the main power cord.

General Instructions

Read Safety Instructions: Unpack the stirrer carefully. Ensure the stirrer is not damaged and verify the contents against the packing list. If the stirrer is damaged or the contents do not correspond with the packing list, notify your supplier immediately.

Read the manual carefully! Make sure that all users read this manual thoroughly before operating the stirrer. For further questions or inquiries, please contact:

Caframo Limited

501273 Grey Road 1
Georgian Bluffs ON N0H 2T0, Canada


Tel: 519-534-1080
Toll Free: 800-567-3556 (within USA and Canada)
Fax: 800-209-6786
Website: www.CaframoLabSolutions.com
E-mail: labsales@caframo.com

Approvals

Independently tested and verified to conform to the following standards:

IEC 61010-1:2001 (2nd Edition)
EN 61010-1:2001 (2nd Edition)
Safety Requirements for Electrical Equipment for Measurement, Control and Laboratory Use – Part 1

Description

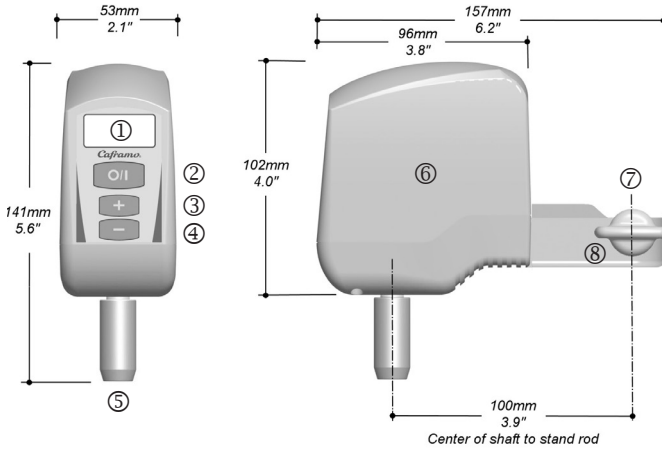- | | |
|--------------------------|-----------------|
| ① Four-digit LED display | ⑤ Collet |
| ② Button - power on/off | ⑥ Housing |
| ③ Button - Increase rpm | ⑦ Clamp opening |
| ④ Button - Decrease rpm | ⑧ Clamp knob |

General Operating Instructions

The stirrer must be securely mounted as specified in the installation instructions. Ensure the mixing shaft and blade are securely attached. The stirrer is now ready for use to mix liquids.

Button Operation Functions

Power Button - Turns the stirrer motor on and off. When initially turned on, stirrer will spin at 50 rpm.

Increase Button - Increases the speed.

Decrease Button - Decreases the speed.

OL Indicates an overload condition. The load is too great for the stirrer and must be reduced. The display will flash "OL" five times and the unit will turn off. To reset stirrer turn power back on.

Technical Data

Model Number	BDC250
Speed range	50 to 2500 rpm
Maximum torque	10 Ncm
Mechanical output power	1/100 hp, 7 watts
Input to power supply	AC 100 to 240 volts, 50/60 Hz (must use supplied power cord)
Electrical output from power supply to Stirrer	DC 24 volts, 0.5 amps
Weight - Stirrer only	0.37 kgs (0.8 lbs)
Weight - Total package	1.47 kgs (3.2 lbs)
Maximum volume	2 litres, water-like viscosity
Molded housing	Chemical-resistant, high-impact nylon. Recyclable (PA)
Keypad	Waterproof, chemical-resistant polyester
Power cord length	1.8 m (6 ft)
Stainless steel collet	Fits 6 mm shaft
Led display accuracy	± 10 rpm or $\pm 3\%$, whichever is greater, 4 digit

Environmental Operating Conditions

The BDC250 stirrer must operate in the following conditions:

Indoors

1. Altitudes up to 2000 m (6500 ft.).
2. Temperatures from 5°C to 40°C (41°F to 104°F).
3. Maximum relative humidity 80% for temperatures up to 31°C (88°F) decreasing linearly to 50% relative humidity at (104°F).
4. Pollution degree 2 in accordance with IEC 664.
“Normally only non conductive pollution occurs. Occasionally, however, temporary conductivity caused by condensation must be expected.”

Installation in the Laboratory (BDC250MU1 includes motor only)

Stand Assembly

Remove lock nut. Insert the threaded end of the stand rod into the stand base. Turn rod clockwise to tighten. Secure with lock nut.

Stirrer Installation

Slide the clamp opening over the top of the stand rod. Lower to desired height. Turn clamp knob to securely tighten the stirrer onto the stand rod. Connect supplied power cord to stirrer. Connect power cord to a grounded electrical outlet.

Impeller Assembly

Add impeller blade(s) to impeller shaft. Tighten screw to secure.

Impeller Installation

Insert impeller shaft into collet opening until it stops. Tighten set screw.

Cleaning and Service

Cleaning

Using a mild soap and water solution, wipe the housing and control panel of the stirrer with a damp cloth. Do not use chlorine bleach, chlorine-based cleanser, abrasives, ammonia, steel wool, or scouring pads with metal content or similar harsh solvents or abrasives. These may damage the surface of the stirrer.

Service

Do not attempt to service or repair stirrer. If the stirrer housing is opened, the warranty becomes void. Contact Supplier for return authorization and return instructions. Clean the stirrer so that it is free from any materials which may be harmful. Provide an SDS or other safety data sheets for all chemicals used in conjunction with this stirrer where appropriate. Place the stirrer and its parts into the original packing to protect and prevent damage during transport. Follow shipping directions provided by Caframo when return is authorized and RMA issued.

CAUTION:

 Failure to clean, service, and handle the stirrer as outlined above can cause it to fail which could impair the safety of the user.

Warranty and Product Liability

Warranty

Caframo warrants the herein described product of its manufacture for three years from receipt of the equipment by the purchaser, against defects in material and workmanship. Warranty registration is available on the Service and Support page of our website (www.CaframoLabSolutions.com) or by completing the supplied warranty registration card. Mail or fax back as indicated.

This limited warranty covers parts and labour. If the stirrer housing is opened, the warranty becomes void.

This warranty may only be altered by a specifically published amendment. No individual has authorization to alter the provisions of this warranty policy or its amendments. Expendable items are not covered by this warranty. Damages due to corrosion or accidental breakage are not covered.

Product Liability

Under no circumstances shall Caframo be liable for indirect, consequential, or special damages of any kind.

Caframo reserves the right to make technical changes without prior notice.

Accessories

The following accessories are available for purchase.

Shaft and Impellers - Part No. S100

- (1) Mixing shaft, 186 x 6.35 mm
- (1) Radial blade, 12.7 x 30.5 mm
- (1) Axial blade, 12.7 x 30.5 mm

Stand - Part No. A215

- (1) Base
- (1) Hollow stand rod with lock nut, 495 x 9.5 mm

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